

HOW TO USE CANON GEL CAMERA

Gel Set-up

- (1) Get gel in place.
- (2) Take off gloves and then position the camera.

Please DO NOT get Ethidium Br all over the camera and the keyboard and mouse. Thanks!

Camera Set-up

- (1) Make sure camera has power.
- (2) Make sure little end of USB is plugged into camera in proper port, and the large end is plugged into the computer's USB (could be the monitor which has a USB bus.)
- (3) Turn camera on.

Taking pictures

- (1) Start Adobe Photoshop.
- (2) Menu bar: FILE→IMPORT→"WIA Canon PowerShot S3 IS..."
- (3) A little window pops up where you select the take picture option.
- (4) Import pictures into Photoshop to crop, contrast adjust, etc.
- (5) Save Image to your folder or upload to your SFTP account.
- (6) Delete the images on the camera memory card using windows explorer. I.e., goto computer icon right click on "Canon PowerShot S3 IS"; select "Explore"; and then find images and delete them.

Clean-up

- (1) Remove camera from gel box with gloves off.
- (2) Put gloves back on.
- (3) Remove gel; clean up and wipe down the light box.